

TYRODE (1 Liter)

NaCl	8 g	137 mM
KCl	0.2 g	2.68 mM
26.5% CaCl ₂ · 2H ₂ O	1 mL	1.8 mM
4.42% NaH ₂ PO ₄ · H ₂ O*	1 mL	0.32 mM
Glucose	1 g	5.56 mM
NaHCO ₃	1 g	1.16 mM

Add distilled water up to 1000 mL pH=7.4

* Na₂PO₄ · 2H₂O is 5% *

Filter and **oxygenate** before use.

Keep **refrigerated** between use.

TYRODE (2 Liters)

NaCl	16 g	137 mM
KCl	0.4 g	2.68 mM
26.5% CaCl ₂ · 2H ₂ O	2 mL	1.8 mM
4.42% NaH ₂ PO ₄ · H ₂ O*	2 mL	0.32 mM
Glucose	2 g	5.56 mM
NaHCO ₃	2 g	1.16 mM